


ARC - Assay Ready Cells

before the start:


- Carefully read the protocol.
- Equilibrate all media and buffer to room temperature.
- Cryopreserved cells are very fragile. Keep the cells on dry ice before thawing and process quickly!

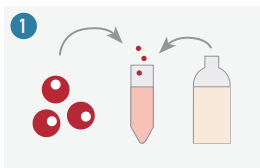

Thaw a vial of Assay Ready Cells in a water bath at **37°C** for **3min**. Transfer the cells into a **15ml** centrifuge tube prepared with **8ml** of Recovery Buffer.



Centrifuge for **3min** at **100xg** and carefully aspirate the supernatant. Resuspend the cell pellet in Assay Medium to your desired cell concentration. Mix gently and let the cells recover for **30min** at room temperature.

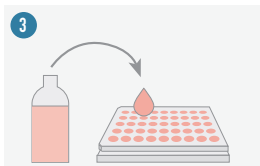

The Cells are **assay ready** now and can be processed according to your assay protocols.

kit content

- ARC - Assay Ready Cells (1x vial)
- Recovery Buffer (1x bottle)
- Assay Medium (1x bottle)

storage

- Store Assay Ready Cells in liquid nitrogen (below -140°C)
- Store all reagents and media at -20°C

limited product warranty

This warranty limits acCELLerate's liability to the replacement of this product. No other warranties of any kind, expressed or implied, including without limitation, implied warranties of merchantability or fitness for a particular purpose, are provided by acCELLerate. acCELLerate shall have no liability for any direct, indirect, consequential, or incidental damages arising out of the use, the results of use, or the inability to use this product. The product is for research only and must not be used for diagnostic or therapeutic purposes.

limited use license

ARC - Assay Ready Cells are sold under a limited use license and shall only be used instantly for an assay application. No further expansion, passaging or refreezing of the cells is permitted. When breaking the sealed bag of the cells the user is explicitly accepting the terms of this limited use license.

Version 1.5

© acCELLerate GmbH (2020)

EUROPEAN OFFICE

+49 (160) 987 577 56

acCELLerate GmbH
Osterfeldstraße 12-14
22529 Hamburg
Germany

please@accelerate.me
www.accelerate.me



US OFFICE

+1 (732) 698 34 04

acCELLerate GmbH
1 Jill Court, Bldg. 16 Unit 10
Hillsborough, NJ 08844
USA